



Orion Clean Card® PRO

Orion Clean Card® PRO

Clean surface? Wipe it and you'll know!

- detects protein residues on surfaces
- simple procedure:
moisten – wipe – interpret
- colour change indicates an unclean surface
- results available in only 30 seconds

HYGIENE

Orion Clean Card® PRO

Food manufacturers and caterers must take every precaution to ensure food safety and good food quality. Adequate cleaning is an essential part of food hygiene practice to prevent food poisoning pathogens and other harmful micro-organisms from spreading between food, surfaces and equipment.

Orion Clean Card® PRO is a new-generation hygiene test for monitoring the effectiveness of cleaning the surfaces. The test detects protein residues, the presence of which indicates inadequate cleaning.

Each Clean Card PRO test device contains a pad impregnated with test reagent. The pad is located under the perforated part of the test device. When proteins are brought into contact with the test device, a visually detectable colour change from yellow to green or blue will occur.

In addition to food safety Clean Card PRO can be used also in other application areas like:

- Monitor the function of your dish washers by checking the cleanliness of washed dishes
- Check the cleanliness of critical control points in hospital wards
- Ensure that your instruments are free of protein residues before autoclaving them

Features

- simple procedure: moisten – wipe – interpret
- colour change indicates an unclean surface
- results available in 30 seconds
- easy to interpret
- long shelf-life
- storage at room temperature
- can be performed by anyone in the staff
- no costly investment
- no instrument needed
- available in two package sizes



Orion
Clean Card® PRO

ORION
DIAGNOSTICA

Orion Diagnostica Oy, P.O. Box 83, FI-02101 Espoo, Finland
Tel. +358 10 4261, Fax +358 10 426 2794, www.oriondiagnostica.com

Orion Clean Card® PRO

Contents of the kit	Cat. No.
<ul style="list-style-type: none">• 25 test devices• Plastic bottle• Spray cap for the bottle• Model chart• Instructions for use	133148
Contents of the kit	Cat. No.
<ul style="list-style-type: none">• 50 test devices• Model chart• Instructions for use	133149
Storage	18–25 °C
Patents	pending

Orion Clean Card PRO is a registered trademark of Orion Diagnostica Oy.

It's so easy to use!

Fill the bottle provided in the kit with clean water. Screw the spray cap on the bottle. Interpretation: If the perforated part remains yellow, the sampled surface is clean. If the colour of the perforated part changes from yellow to green or blue, the sampled surface is not clean. The darker the blue colour becomes, the more protein there is on the sampled surface.

